# MR-VP BROTH

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pH = 7,0 +/- 0,2

## PREPARATION

Suspend 17 gr of powder in 1 litre of distilled or water. Heat until completely dissolved. Dispense in tubes and sterilize in the autoclave at 121°C for 15 minutes.

## USE

MR-VP BROTH is a medium used to differentiate strains of coliform bacteria by the Methyl Red (MR) and Voges Proskauer (VP) tests. *Escherichia coli* and other MR positive bacteria ferment glucose with high acidity production. *Enterobacter aerogenes* and other MR negative bacteria produce less acid reaction and on continued incubation became more alkaline as the acid metabolites are transformed in carbonates and carbon dioxide. The methyl red test is carried out by adding 5 drops of methyl red solution at 0,02% to 5 ml of broth culture in MR-VP MEDIUM incubated at 30°C for 5 days. A positive reaction is indicated by a distinct red color while a negative reaction is indicated by a yellow color. The Vooges-Proskauer reaction differentiate the bacteria fermenting glucose into acids from those which ferment it into acetyl methyl carbinol which gets oxydated into diacetyl by the oxygen present in an alkaline environment following the addition of NaOH. Diacetyl reacts with alpha naphthol developing a pink-red color. The test is carried out by adding the following solutions to 1 ml of broth culture in MR-VP MEDIUM after 24-48 hours of incubation at 37°C:

- 1) 0,2 ml of 40% NaOH watery solution
- 2) 0,6 ml of 5% alpha naphthol alcoholic solution

The positive reaction is indicated by a change in color from pink to red 2-4 hours following the addition of the reagents.

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**BIBLIOGRAPHY**

1. APHA (1963). *Diagnostic Procedures and Reagents*, 4<sup>th</sup> edition.
2. Barrit M.M. (1963). *The Intensification of the Voges Proskauer Reaction by the Addition of Alphanaphthol*. *J. Pathol. Bacterio.*, 42, 441.
3. Edwards P.R. and Ewing V.H. (1955). *Identification of Enterobacteriaceae* Minneapolis: Burgess Publishing Company.