TRYPTIC (CTA) MEDIUM

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PREPARATION

Suspend 28,5 gr of powder in 1 litre of distilled or deionized water. Heat until completely dissolved. Dispense in 10 ml tubes and sterilize in the autoclave at 121°C for 15 minutes. Cool to 50°C. If desired aseptically and a sterilized solution for filtration of the appropriate carbohydrate so as to obtain a final concentration of 0,5-1%.

USE

TRYPTIC (CTA) MEDIUM is a medium used for the growth and preservation of pathogenic microorganisms and to carry out mobility test. By adding carbohydrates the bacteria fermentative metabolism can be observed. The low concentration of the agar (0,25%) reduces the oxygen diffusion and the carbon dioxide release allowing the growth of the anaerobic bacteria. Furthermore, by inoculating a TRYPTIC (CTA) MEDIUM tube by threading, the microorganisms with mobility will show growth from the inoculation line. Fermentation of the added carbohydrates, determines an acidification of the medium and a color change of the red phenol indicator.

CULTURAL CHARACTERISTICS AFTER 24-48 HOURS OF INCUBATION AT 37°C

Microorganism	Growth	Mobility	Acid production
<i>E. coli</i>	good	+	+
<i>N. gonorrhoeae</i>	good	-	+
<i>N. meningitidis</i>	good	-	+
<i>Strept. pneumoniae</i>	good	-	+

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