CLED ANDRADE AGAR

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PREPARATION

Suspend 36,3 gr. of powder in 1 litre of distilled or deionized water. Heat until completely dissolved. Dispense and sterilize in the autoclave at 121°C for 15 minutes.

USE

CLED ANDRADE AGAR is a non selective medium suitable for the estimation of bacteria from urine. Andrade allows a more accurate detection of lactose fermenting bacteria from non fermenting ones. The more or less intense fermentation of the lactose causes a color change in the medium towards different red tones. The unsuccessful fermentation of the lactose causes an alkalization of the medium with consequent color change to green-blue.

CULTURAL CHARACTERISTICS AFTER 18-24 HOURS OF INCUBATION AT 37°C

Microrganismo	Colonies colour
<i>Escherichia coli</i>	red with a pink-red halo
<i>Proteus mirabilis</i>	green-blue with irregular edges
<i>Klebsiella pneumoniae</i>	muculent, grey-green
<i>Staphylococcus aureus</i>	small, yellow
<i>Streptococcus faecalis</i>	small, yellow-orange

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BIBLIOGRAPHY

Mackey J.P. and Sandys G.H. (1965). *Laboratory Diagnosis of Infections of the Urinary Tract in General Practice by Means of a Dip-Inoculum Transport Medium*. Br. Med. J., 2; S 1286.