# LISTERIA UVM 1 BROTH

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## PREPARATION

Suspend 54 gr of powder in 1 litre of distilled or deionized water. Heat until completely dissolved. Dispense in vials and sterilize in the autoclave at 115°C for 15 minutes.

## USE

LISTERIA UVM 1 BROTH is a selective broth for the isolation and cultivation of *Listeria*. The USDA-FSIS procedures recommend LISTERIA UVM 1 BROTH as a pre-enrichment broth for the identification of *Listeria* in meat. The selective agents contained are acriflavine and nalidixic acid.

Procedure:

- 1) Homogenize 25 gr of specimen in 225 ml of LISTERIA UVM 1 BROTH and incubate at 30°C for 24 hours;
- 2) Transfer 0,1 ml of UVM 1 BROTH broth culture in 10 ml of Listeria UVM 2 Broth and incubate at 30°C for 18-24 hours;
- 3) Take a loopful of broth culture from Listeria UVM 2 Broth and streak selective and differential medium for *Listeria* (*Listeria Aesculin Agar*, *Listeria Palcam Agar*) placed on Petri dish. Incubate at 30°C or 37°C for 24-48 hours;
- 4) The suspect colonies must be submitted to Gram coloring, catalysis test, mobility examination and identification biochemical tests.

## CULTURAL CHARACTERISTICS AFTER 24 HOURS OF INCUBATION AT 30°C

Microorganism	Growth
<i>Listeria monocytogenes</i>	good to excellent
<i>Staphylococcus aureus</i>	partially to completely inhibited
<i>Escherichia coli</i>	partially to completely inhibited

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